PCRFast[®] Mandel

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09/07/2008

PCRFast[®] Almond

Realtime (SYBR[®] Green) and gel detection

PCR - test for the detection of almond

Type No.: IF / AG1002

In-vitro test

Storage: 2 - 8 °C (35.6 - 46.4.°F)

Brief information

Easy-to-use molecular biological test (PCR) for the detection of almond (*prunus dulcis*, detection of the lipid-transfer-protein I - gene with 98bp) in food and pharmaceutical products. The test kit contains 96 reaction vials. All reaction vials contain a specific primer pair, and in addition 48 reaction vials (red-coloured) contain specific almond DNA for the PCR positive control and for the inhibition control.

Usage

For the general usage of the kit see "PCRFast[®] - General Information". Please read the instruction carefully.

Cycler profile

10 min	95 °C (203 °F)	45 cycles
15 sec	95 °C (203 °F)	
60 sec	61 °C (145.4 °F)	

The sequential identity of the amplified DNA (98bp) could be verified by a restriction analysis with *Fnu4H*. This results in two fragments with a length of 39bp and 59bp.

Sensitivity

The limit of detection is < 10 copies.

Specificity

PCRFast[®] Almond is specific for almond and nectarine. The following species have been tested with 100 ng DNA for cross reactivity:

species		species		species		species	
almond	+	nectarine	+	sesame	-	pine	-
cabbage	-	orange	-	celery	-	dill	-
anise	-	lupine	-	pea	-	cashew	-
apricot	-	hazelnut	-	walnut	-	pecan	-
macadamia	-	brazil nut	-	pistachio	-	peanut	-
corn	-	soy	-	sunflower	-	pumpkin	-
lentil	-	rye	-	buckwheat	-	barley	-
sorghum	-	5-cereal-mix	-	oat	-		

Tab. 1: Specificity PCRFast[®] Almond

+ : band with 98bp

- : no band with 98bp

Lit: Sambrook J, Fritsch E F, Maniatis T

„Molecular cloning. A laboratory manual“, (3rded.), New York: Cold Spring Harbor Laboratory Press.

Certificate of Quality

PCRFast[®] Almond,
complies specification.

Lot: TMA 11111
quality assurance: A.Schuhr

09/07/2008