PCRFast[®] VTEC

Realtime (Sonde)

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PCRFast[®] VTEC

Realtime (Probe)

PCR - test for the detection of shigatoxin producing Escherichia coli (VTEC stx1 and stx2)

Type No.: IF / MR1004

In-vitro test

Storage: 2 - 8 °C (35.6 - 46.4 °F)

Brief information

Simple, molecular biological test (PCR) for the qualitative detection of shigatoxin producing Escherichia coli in foodstuff by gel detection (detection of the genes for stx1 and stx2, probe with Fam reporter and non fluorescence quencher). The test kit has 96 reaction vials, each reaction vial contains specific primers, probes and an internal amplification control (probe with HEX reporter and non fluorescence quencher) for investigating possible inhibiting effects (ITC). 8 reaction vials (red marking) contain additionally VTEC DNA for the PCR positive controls.

Usage

For the general usage of the kit see "PCRFast[®] - VTEC (probe)". Please read the instruction carefully.

Cycler profile

10 min	95 °C (203 °F)	
15 sec	95 °C (203 °F)	
60 sec	60 °C (140 °F)	45 cycles

Sensitivity

The limit of detection is < 10 copies.

Specificity

PCRFast[®] VTEC is 100 % specific for VTEC stx 1 and stx 2. The following species have been tested with at least 2500 copies for cross reactivity:

Species		Species		Species	
VTEC stx1	+	VTEC stx2	+	Escherichia coli	-
Yersinia enterocolitica	-	Legionella pneumophila	-	Campylobacter jejuni	-
Staphylococcus aureus	-	Bacillus cereus	-	Salmonellae	-
Clostridium perfringens	-	Listeria monocytogenes	-	Legionella erythra	-

Tab. 1: Specificity PCRFast[®] VTEC

+ : amplification

- : no amplification

ISO 22174

Microbiology of food and animal feeding stuffs - Polymerase chain reaction (PCR) for the detection of food - borne pathogens - General requirements and definitions

Certificate of Quality

PCRFast[®] VTEC,
complies specification.

Lot: **RECD 11111**
quality assurance: A.Schuhr

20/10/2008